Zeiss 880 Training Notes Zen 2.3





Logging into the 880

- Log into the computer with your UCHC or CAM credentials; be sure you are logging into the proper domain.
- Double click on the Zen icon, the one labeled with "AUT-Zeiss LSM 880" or similar text.

The idea elan tracking and billing software will initialize after clicking on the Zen icon.

 Enter your UCHC email. Once it has been recognized, it will bypass the password requirement and proceed to the session confirmation window.





Starting Zen with idea elan

 Select your scheduled appointment and/or click the login button as directed by the software.



• Confirm your appointment and the Zen software will be initialized.



Starting Zen 2.3

Select **Start System** from the Zen 2012 login window. This will start up the hardware and software.

Image Processing is for using the system in an offline mode and this will ignore all the hardware components.

	ZEN	
1000		
Boot Status		
Start System		Image Processing
Start System		Image Processing

Introducing the Zen Workspace

The Zen workspace is designed in a tabular fashion allowing you to step you through initially looking at your specimen via the Oculars tab, in the control panel, to acquiring images via the Acquire tab. The remaining tabs add offline Processing and Maintenance options.



Panels can be moved around `to customize the workspace.

Introducing the Zen workspace

Your imaging configuration and your work space configuration can be stored separately.

In addition, your imaging configuration and acquisition parameters can easily be retrieved from the **Reuse** function when you have a previously saved image loaded in the workspace.





Getting Light to the Oculars

Locate Tab

- 1. Click on the objective icon to access the list of possible objectives. Be sure to note the immersion type; air, oil or water. Please do not change from oil to air/dry or from water to oil and oil to water without cleaning your specimen first.
- Select your light source from: Transmitted ON Transmitted OFF DAPI FITC

Rhodamine

3. Be sure to turn the Reflected Light "Off" when you are finished looking at your specimen.



Getting all the tools

Acquisition Tab

In order to view all the options available via the software, be sure to go to:

- 1. View>Show All (global)
- 2. Select >Show all Tools.

Make sure to do this each time you use the software otherwise the required features, like the lasers, will not be accessible to you.

Open/close the submenu with the arrow.

Click the arrow to detach/attach panel.



Acquisition Tab

1. Experiment Manager

The manager allows you to save your scan configuration settings for your experiment and reload the settings for a quick start. This method does not record the acquisition parameters.

- 2. Scan Control Tool Bar
 - AF (Find Focus) computer determines focus based on contrast settings.
 - Set Exposure the computer automatically calculates the gain and black level settings.
 - Live initiates a fast scan regardless of the scan speed and continuously scans until Stop is selected.
 - Continuous continuously scans according to the selected scan speed until Stop is selected.
 - Snap performs a single scan according to the selected scan speed.



Using Smart Setup

Smart Setup

The best and easiest way to get started is to use the Smart Setup tool.

- 1. Click on the down arrow to open the dye selection menu.
- 2. Select your dye from the drop down menu. Repeat this procedure to select additional dyes.

Depending on your imaging needs, the Smart Setup will offer you up to four different configurations for imaging your sample:

- Fastest
- Best Signal
- Best Compromise
- Linear Unmixing.

Whatever configurations you are presented with, you can always manually modify the configuration in the Light Path panel.



Smart Setup - Fastest

Acquires the image with the simultaneous excitation and detection of one or multiple probes. This option produces a single image, with a single scan however it may result in bleed through between the different channels.



Smart Setup – Best Signal

Acquires the image with the sequential excitation and detection of the different probes. Each probe is scanned separately and the final image is the result of an overlay of the individual channels. This option avoids bleed through between the different channels however multiple scans are require to produce the image which can lead to bleaching or other laser induced damage.



Smart Setup – Smartest (Line)

- Smartest acquires multiple channels at the same time while quickly toggling laser lines to acquire a second track.
- There are limitations to this option, especially if you are acquiring 4 different probes.
 - a) You must you use the same dichroics between the two tracks.
 - b) You can't send two probes, of different intensities to the ChS detector.



Smart Setup – Linear Unmixing

This option creates a confirmation for computationally subtracting bleed through between channels. This can be performed after you acquire individual, pure spectral profiles of the probes you are using.



Using the Advanced Scanning Modules

Advanced scanning modules are available as a single option or they can be used in combination with multiple modules. The modules are initiated via the Start Experiment button.

- 1. Time Series collect a series of images based on a specific time interval and duration.
- 2. Z-Stack collect a series of optical sections that can be reconstructed as a 3D image.

cxperimenting					u Multidimonsional /	Acquicition		
4 color 4 dete	ctors			- H 🖬 🗇	• Waldamensional #	Acquisition		
Y Smort Sot				(🔽 🗏 Z-Stack		√ s	how All [
7 aniai caec	up		snow all loois	New	First/Last			
AF	0	_		Ó		Cotta	ct 10.00	
Find Focus	et Exposure	Live	Continuous	Snap		Set La	50 10.00	
7-Stack	21 Slires						20.00	ım
Time Series	21 Jiices						21	\$
Bleaching							1.00	13
Tile Scan								i lim
Positions			Start Exp	eriment				μm
_ negions							Interval	Slice
				-	-14.1	3 Set Fil	rst -10.00	:
🕨 🦞 Laser					Position (um) 0.00			
F Light Pa	th		🗸 Sho	WAII	rosition (pm) - 0.00			



 • • • Time Series 	🗸 Show All 🛛 📝
Cycles -[]	10 🛟
	<u>10</u>
G Interval Time	
Interval Time not defined	
Marker	
Marker not defined	📄 🗄 🖻
+ -	
🗢 Start	
Mode Manual 👻	📄 Pre-Scan
Trigger Out None 🔫	
• End	
Mode Manual 👻	
Trigger Out None	
Pause	

Using the Advanced Scanning Modules (cont.)



Bleaching – module for FRAP (fluorescent recovery after photobleaching.)

	Show all 🕜
Bleach settings not defined	
Start Bleaching after # scans	
Repeat Bleach after # scans	
Iterations 1	•
Different scan speed	
Safe bleach for GaAsP	
Different Z Position [µm]	
Trigger in None 🔻 Trigger	out None 🔻
Zoom Bleach (fast, less accurate)	
Excitation of Bleach	
Use different Settings for different RO	
ROI# all	
R01# all	-
ROI# all 405 458 488 514 561 633	
R01# all 405 458 488 514 561 633 Test Bleach	



Tile Scan – create larger images by tiling/stitching smaller images together.







Using the lasers

The lasers that are available for use are:

Diode 405 nm Argon 458,488,514 nm DPSS (Diode-pumped solid state) 561 nm HeNe (Helium-Neon) 633 nm

The Argon, has a start-up and shut-down procedure. This can be monitored by opening the **Laser Properties** panel.

The red highlighted color indicates the laser is not available for use. (Please note this color will go away, it will not turn green, when the laser is warmed up.)

👻 🂡 Laser			
Laser	Laser Lines [nm]]	Power
Argon	458, 488, 514		(On 🔹
🔺 Diode 405-30	405		
📤 DPSS 561-10	561		On 🔻
📤 HeNe633	633		On 🔻
Laser Properties			
Wavelength	458, 488, 514 nm		
Status	Busy		
Tube Current	4.4 A		
Power	Optimal 🔻	32 %	

Looking at the Light Path

1. Tracks

- Create and/or delete tracks which contain the scanning configuration for each probe or combination of probes that are scanned. The configuration consists of the laser line, excitation dichroic, detector and emission range that is captured. If you use the Smart Setup application, your tracks will automatically be configured for you.
- 2. Switch track every
 - Line tracks are switched line by line. Laser line, laser intensity and channels can be changed between tracks.
 - Frame tracks are switched frame by frame. Laser line and intensity, filters, beam splitters, gain and offset, and pinhole setting can all be changed between tracks.
 - Frame Fast scanning procedure can be made faster. Only the laser line intensity and Amplifier offset can be changed between tracks.



Imaging Setup and Light Path

- Track defines the fluorophore(s) that is(are) being detected. A maximum of four tracks can be created which specify the dye, laser excitation line and the emission detection range.
- Visual representation of excitation line(s) and the emission profile(s) which you can set by adjusting the slider bars or alternatively you can enter the range at the drop down arrow.



Imaging Setup and Light Path

1. Tracks are defined by dye, pseudo color, detector, and emission range. The check mark indicates the track is enabled and will scan upon activation.

> Detectors – select from standard PMTs, Ch1 and Ch2 or the QUASAR detector ChS 1-8, and the transmitted light detector ChD (not confocal).

 Select laser wavelengths, either visible or invisible (UV), and their attenuation values (% transmission) and the main dichroic beam splitters.



Acquisition Mode

- 1. Objective change objectives, this is an alternative to going to the locate tab.
- 2. Scan Mode, Frame Size and Line Step
 - Scan Mode Frame, Line, Spot
 - Frame Size defaults to 512 x 512. Select Optimal for appropriate number of pixels depending on numerical aperture and excitation wavelength.
 - Line Step scan every line or less than every line.
- 3. Speed adjust scan speed. A higher speed with averaging yields results with the best signal-tonoise ratio. Use slower scan speeds for better images, but be aware of photobleaching.



Acquisition Mode (cont.)

- Averaging improves image by increasing the signal-to-noise ratio. Select from Line or Frame averaging and select the number to average by. Averaging helps to reduce photobleaching.
- Bit Depth select from 8 bit (256 gray levels), 12 bit (4096 gray levels), or 16 bit (65536 gray levels). Be sure your software can handle the information. 12 and 16 bit images are recommended for publication and quantitative measurements.

Direction – x-y scan direction or bi-directional

6. Scan Area – set optical zoom setting and /or rotate laser orientation.



Working with Channels

The channels dialog allows you to select an individual track and adjust the corresponding settings to improve the image quality.

- Tracks as defined from the light path. Be sure to highlight the channel you wish to modify. If a track has a "check mark", it is active and it will be scanned.
- 2. Lasers enable or disable and modify laser power.
- Pinhole set pinhole to 1 AU (airy unit) to achieve best z resolution based on the excitation wavelength and objective numerical aperture. Be sure to match the section thickness when scanning multiple channels.



Working with Channels(cont.)

- 4. Gain, Digital Offset, Digital Gain
 - Gain (Master) amplifies the voltage on the photomultiplier tube; regulates brightness of the image.
 - Digital Offset sets background/threshold level or dark current of the image.
 - Digital Gain post acquisition enhancement of the gain signal.

Tracks Channels ✓ Track 1 A488 ✓ Track 2 DAPI ✓ A568 ✓ ✓ Track 3 A647 ✓ ✓ ★ ✓ Track 3 A647 ✓ ✓ ★ ✓ Track 3 A647 ✓ ✓ ★ ✓ A ● Track Configuration not defined ✓ Track 2 - LSM Lasers ✓ ● Lasers ✓ ● ● ● 405 nm 2.0 ● ● ● ✓ ✓ ● ● ● ● S61 nm 2.0 ● ● ● ● Jog Airy Units ± 1.9 µm section IAU max ● ● ● Digital Offset ● ● ● ● ● Digital Offset ● ● ● ● ● Digital Offset ● ● ● ● ● ●	🗢 📇 Ch	annels			~	/ Show all	B
Tracks Channels ✓ Track 1 A488 ✓ Track 2 DAPI A568 ✓ ✓ Track3 A647 ✓ Track4 A647 ✓ Track3 A647 ✓ Track4 A647 ✓ Track5 A647 ✓ Track 2 Lasers ✓ A + Expand All Collapse All Track 2 - LSM Lasers Lasers ✓ ✓ 551 nm 2.0 ✓ 561 nm 2.0 Pinhole 34.1 1.39 Airy Units ≜ 1.9 µm section 1 AU max DAPI Gain (Master) 435 Digital Offset 0 2 Digital Offset 3 2 Digital Offset 3 2 Digital Offset 3 2 Digital Gain 1.0 2							
✓ Track 1 A488 ✓ Track 2 DAPI A568 ✓ Track3 A647 ✓ Track 2 Losers ✓ A + 10 Expand All Collapse All Track 2 - LSM Lasers Lasers ✓ 458 ✓ 405 458 ✓ 561 nm 2.0 ° ✓ 561 nm 2.0 ° ✓ 561 nm 2.0 ° ✓ 1.39 Airy Units ≤ 1.9 µm section 1.4U max DAPI Gain (Master) 1.33 ° Øigital Offset 0 0 Øigital Offset 3. ° 1.0 ° Øigital Offset 3. ° 1.0 ° Øigital Gain 1.0 ° 1.0 °	Trac						
✓ Track 2 DAPI A568 ✓ Track3 A647 ✓ Track3 A647 ✓ Track3 A647 ✓ Track Configuration not defined Track Configuration not defined Track 2 - LSM Lasers ✓ 405 458 405 458 405 458 405 458 405 435 561 nm 2.0 Pinhole 34.1 1.39 Airy Units ≜ 1.9 µm section DAPI Gain (Master) 553 Digital Offset 0 2.0 Digital Offset 3 2 Digital Offset 3 2 Digital Offset 3 2 Digital Offset 3 2 Digital Gain 1.0 2							
A568 ✓ Track3 A647 ✓ ↑ + ∅ Expand All Collapse All Track Configuration not defined Track 2 - LSM Lasers 405 405 458 405 458 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 405 435 561 1.9 µm section 1.39 Airy Units ≜ 1.9 µm section DAPI Gain (Master) Digital Offset 0 Digital Offset 0 Digital Offset 3 Digital Offset 3 Digital Gain 1.0						-	
✓ Track3 A647 ✓ + Expand All Collapse All Track Configuration not defined ✓ H Track 2 - LSM ✓ ✓ Lasers ✓ ✓ ✓ 405 488 514 561 633 405 A647 ✓ ✓ ✓ Pinhole ✓ ✓ ✓ ✓ 1.39 Airy Units ≜ 1.9 µm section 1.40 max DAPI Gain (Master) ✓ ✓ Digital Offset 0 ✓ ✓ Digital Offset 0 ✓ ✓ Digital Offset 0 ✓ ✓ Digital Gain ✓ ✓ ✓						-	
Image: space of the							
Track Configuration not defined Image: Configuration not defined Track 2 - LSM Lasers Image: Configuration not defined 405 458 405 458 405 458 405 458 405 405 405 488 561 633 405 435 1.39 Airy Units ≤ 1.9 µm section AAF68 Gain (Master) 0 1.0 A568 Gain (Master) 0 1.0 0 1.0							
Track 2 - LSM Lasers 405 458 488 514 561 633 405 nm 2.0 2.0 2.0 2.0 2.0 ▲ 561 nm 2.0 34.1 2.0 2.0 2.0 Pinhole 34.1	Track Co					ដា 🕯	
Track 2 - LSM Lasers ✓ 405 488 514 561 405 nm 2.0 ° ▲ 561 nm 2.0 ° Pinhole 0 34.1 ° 1.39 Airy Units ≜ 1.9 µm section 1.4U max DAPI Gain (Master) 0 553 ° Digital Offset 0 0 20 Digital Offset 0 553 ° 1.0 ° Digital Offset 0 3 20 Digital Offset 0 1.0 ° 3 20							
Lasers ↓ 405 458 488 514 561 633 405 nm 1 2.0 2 ▲ 561 nm 2.0 2 Pinhole 34.1 2 1.39 Airy Units ≜ 1.9 µm section 1AU max DAPI Gain (Master) 553 2 Digital Offset 0 2 Digital Gain 1 1.0 2 A568 Gain (Master) 435 2 Digital Offset 0 3 2 D	Track						
Lasers 405 458 488 514 561 633 405 nm 2.0 2.0 2.0 2.0 ▲ 561 nm 2.0 2.0 2.0 2.0 ▲ 561 nm 2.0							
405 458 488 514 561 633 405 nm 2.0 ° ▲ 561 nm 2.0 ° Pinhole 0 34.1 ° 1.39 Airy Units ▲ 1.9 µm section 1 AU max DAPI Gain (Master) 0 553 ° Digital Offset 0 ° Digital Gain 1.0 ° 435 ° Digital Offset 3 ° ° Digital Offset 0 1.0 °	Lasers						
405 nm 2.0 : 561 nm 2.0 : Pinhole 1.39 Airy Units ≜ 19 µm section 1 AU max DAPI Gain (Master) Digital Offset Digital Gain 1 0 : A568 Gain (Master) Digital Offset Digital Offs							
▲ 561 nm 2.0 ‡ Pinhole 34.1 ‡ 1.39 Airy Units ▲ 1.9 µm section 1 AU max DAPI Gain (Master) 1 553 ‡ Digital Offset 0 ‡ Digital Gain 1.0 ‡ A568 Gain (Master) 435 ‡ Digital Offset 3 ‡ Digital Offset 1.0 ‡ Digital Offset 1.0 ‡	40						
Pinhole 34.1 ÷ 1.39 Airy Units ≜ 1.9 µm section 1 AU max DAPI Gain (Master) 553 ÷ Digital Offset 0 ÷ Digital Gain 1 1.0 ÷ A568 Gain (Master) 435 ÷ Digital Offset 3 ÷ Digital Gain 1 1.0 ÷	A 56						
1.39 Airy Units ▲ 1.9 µm section 1 AU max DAPI Gain (Master) ● 553 ÷ Digital Offset ● 0 ÷ Digital Gain ● 1.0 ÷ A568 Gain (Master) ● 435 ÷ Digital Offset ● ● 3 ÷ Digital Offset ● ● 1.0 ÷	Pinhole						
1.39 Arry Units 1.9 µm section 1 AU max DAPI Gain (Master) 1 553 1 Digital Offset 0 1 1 1 Digital Gain 1 1 1 1 A568 Gain (Master) 435 1 Digital Offset 1 3 1 Digital Gain 1 1 1	11111010						
DAPI Gain (Master) 553 \$ Digital Offset 0 \$ Digital Gain 1.0 \$ A568 Gain (Master) Digital Offset 435 Digital Offset 3 Digital Gain 1.0	1.39	Airy Units ≙	1.9 μm se	tion	Ŀ	LAU ma	эх
A568 Gain (Master) - 0 : Digital Gain - 1.0 : Digital Offset 3 : Digital Gain - 1.0 :	DAPI	Gain (Maste	er) —	-1		53	÷,
A568 Gain (Master) 1.0 Digital Gain - 435 Digital Offset		Digital Offs	et —				
A568 Gain (Master) 435 Digital Offset			- 0			0	-
A568 Gain (Master) 435 Digital Offset		Digital Gali	n -ll			.0	•
Digital Offset	A568	Gain (Mast	er) —	-0	- 4	35	÷
Digital Gain -() 1.0 🗘		Digital Offs	et		-1-3		
					- 1	<u>^</u>	•
		Digital Gair	n -1)		1	.0	•

Using the Range Indicator

- Remember, when acquiring your images, you want to be within the dynamic range of the detector. This is especially critical when collecting images that will be used for quantitative analysis.
- Use the Range Indicator to adjust the gain and offset. Blue= minimum, Red= maximum (0 and 255 for 8 bit image, 0 and 4095 for 12 image, 0 and 65536 for 16 bit image.)
- Adjust Detector Gain and Amplifier Offset to remove red and blue pixels. Minimum should be ~ 10-20 gray levels above 0 and Maximum should be 10-20 gray levels below 255 for an 8 bit image. You can scale this accordingly for 12 and 16 bit images.

Ch2



Using the Focus Tool

- The 780 offers the ability to manually control the focus or to use the software for more precise control of the focus.
- Step size set a fixed step size, in microns. Move up or down in the focal plane based on that step size.
- 2. Z-position specifies location of focal plane, with regard to a user set zero position.
- 3. Home zeros the focal position.
- 4. Work returns the objective to the desired focus, set with the previous sample.
- 5. Load lowers the objective so the user can change the sample.



Using the Stage Tool

- The 780 allows you to manually control the stage via the joy stick or to use the software to control the stage.
- 1. x-y coordinates indicates stage position.
- 2. Step move specified step size, in microns.
- 3. Set zero establish a zero point (0,0) on the stage
- 4. Move to zero move to the zero point on the stage.
- 5. Marks generate a list of stored positions, with regards to x, y, and z.



Acquiring a Z-stack



Used to collect multiple images, of different focal planes, within a single file for 3-D analysis/display.

- 1. You can enable the z-stack dialog from advanced features menu.
- 2. Scroll down to access and then open the Z-Stack dialog.

Z-stack Dialog

E 2-states	
First/Last	Center
153	Set Last 11.00 ‡
	Range 20.00 µm
	Slices 21
	Interval 1.00
	Smallest 0.69 µm
	Keep 💿 Interval 🏐 Slice
-13.3	Set First -9.00
Position (um) 0.00	
14 μm 14 μm 14 μ	
Match Pinhole Optimal	Undo X.YZ=1
Match Pinhole Optimal Correction Refractive Index Correction	Undo X.Y.Z = 1
Match Pinhole Optimal Correction Refractive Index Correction Auto Z Brightness Correction	Undo X.Y.Z = 1 1.00
Match Pinhole Optimal Correction Refractive Index Correction Auto Z Brightness Correction Use Correction	Undo X.Y.Z = 1 1.00
Match Pinhole Optimal Correction Refractive Index Correction Auto Z Brightness Correction Use Correction Add	Undo X:Y:Z = 1 1.00 • Load Save
Match Pinhole Optimal Correction Refractive Index Correction Auto 2 Brightness Correction Use Correction Add Remove	Load Save Spline interpolation

- 1. First/Last or Center
 - Establish the stack acquisition parameters.
- 2. Optimize Sectioning and Step
 - Ensure the pinholes are set to the same section thickness for multiple wavelengths.
- 3. Correction
 - Options for adjusting refractive index and/or auto brightness to ensure even illumination through the stack.

2

Defining First/Last



Set the first and last starting positions for the stack and acquire the slices within that range.

- 1. Set Last marks z position furthest away from the specimen, as mounted on the cover glass.
- Range total thickness of stack, in microns.
- Slices number of slices making up the stack
- Interval distance between each slice, in microns
- Smallest calculates the slice to be ½ the zresolution for proper Nyquist sampling.
- Keep Interval/Slice selection if the stack should be based on the interval thickness or a specified number of slices.
- 2. Set First marks z position closest to the specimen, as mounted on the cover glass.
- 3. Position Current z position

Optimize Sectioning and Step



- Match Pinhole for multiple wavelengths, ensures that the pinholes are yielding the same section thickness.
- Optimal calculates each slice to be ½ the z-resolution for proper Nyquist sampling.
- 3. Undo clears selection
- 4. X:Y:Z=1 sets the z interval so the voxel is the same for x, y, and z dimensions.

Z-Stack Display Options

- 1. Ortho orthogonal display, xy, yz and xz options
- 2. Cut display a particular slice of stack and have options for manipulating that slice
- 3. 2.5D layered depth display
- 4. 3D reconstruction of z-stack









3

4

Image Display Options

- 1. 2D x y view of image
- 2. Gallery display of z-stack or time series images.
- 3. Split view multi-channel images as individual and overlay displays.







Advanced Image Display Options

- 1. Profile display pixel intensities along defined regions or lines.
- 2. Co-localization –display spatial overlap of fluorescent regions in different channels.
- 3. Histogram display of image intensities and their frequencies within the image.
- 4. Information summary of image file.





Image Display Features



Dimensions Tab

- 1. Z position step through optical z-stack slices
- 2. Zoom digital zoom feature
- Channels range indicator option, pseudo-color selection, turn on/off channels that are displayed
- 4. Ruse load image acquisition parameters for image being displayed.
- 5. Crop isolates a smaller region of the image and optically zooms the image.
- Positions when you click on the image, a crosshair will be displayed, this will correspond to and generate a list of positions within the stage tool.
- Stage cursor appears as a crosshair in a rectangle. As you click and change position, this will update the center position of the stage for the next image.

Image Display Features



Display Tab

- 1. Channel selection– modify display settings for all or individual channels.
- Min/Max, Best Fit methods for adjusting contrast settings.
- 3. Black/White, Gamma set max and minimum levels and non-linear histogram adjustment.

Image Display Features

- 1. Player controller for reviewing z-stack or time series data.
- 2. Graphics image overlays include scale bars, text and other various drawing elements.







Shut Down Sequence Verified Next User Coming

